

Register your instrument! www.eppendorf.com/myeppendorf



CryoCube® F740h

Operating manual

Copyright© 2020 Eppendorf AG, Germany. All rights reserved, including graphics and images. No part of this publication may be reproduced without the prior permission of the copyright owner.

Eppendorf® and the Eppendorf Brand Design are registered trademarks of Eppendorf AG, Germany.

CryoCube® is a registered trademark of Eppendorf AG, Germany.

Registered trademarks and protected trademarks are not marked in all cases with ® or TM in this manual.

U.S. Patents are listed on www.eppendorf.com/ip

Table of contents

1	Oper	ating ins	structions	7
	1.1	Using t	his manual	7
	1.2	Danger	symbols and danger levels	7
		1.2.1	Danger symbols	7
		1.2.2	Danger levels	7
	1.3	Symbol	ls used	8
	1.4	Version	n overview	8
2	Safet	y		9
	2.1	Intende	ed use	9
	2.2	Warnin	gs for intended use	9
		2.2.1	Devices with flammable refrigerant	10
	2.3	User pr	rofile	10
	2.4	Persona	al protective equipment	11
	2.5	Informa	ation on product liability	11
	2.6	Mainte	nance and repairs	11
	2.7	Electro	magnetic compatibility	12
		2.7.1	Europe	12
		2.7.2	U.S.A	12
	2.8	Warnin	g symbols on the device	13
3			ription	
	3.1	Product	t overview	19
		3.1.1	Front view	
		3.1.2	Side view	20
		3.1.3	Internal view	21
		3.1.4	Interfaces	22
	3.2		es	
	3.3		i	
	3.4			
	3.5	Deliver	y package	
		3.5.1	Device and accessories	
		3.5.2	Documents	
	3.6	Accesso	ories	
		3.6.1	Back-up systems	
		3.6.2	Chart recorder	25
		3.6.3	Racks for ULT upright freezers	26
		3.6.4	Cardboard boxes and box dividers	26
		3.6.5	Eppendorf Storage Box	27
4				
	4.1	3		
	4.2		ng installation	
		4.2.1	Unpacking the device	
		4.2.2	Checking the delivery	
		4.2.3	Transporting the device to the location	
		4.2.4	Setting up the device	31

	4.3 Connecting the device to the voltage supply			32
	4.4	Connec	cting the device to external systems	32
		4.4.1	Remote alarm interface	32
		4.4.2	RS-485 interface	33
	4.5	Switchi	ing the device on	33
		4.5.1	Enabling the back-up circuit	34
		4.5.2	Switching the device on at the mains/power switch	34
	4.6	Basic d	device settings	35
_	_			a=
5	•			
	5.1	•	ng the outer door	
	5.2		g the device	
	5.3		g the outer door	
	5.4		re compensation	
	5.5		ing off the device	
		5.5.1	Disabling the back-up circuit	
		5.5.2	Disconnecting the device from the voltage supply	40
6	Softv	vare		41
	6.1	Overvie	ew of the control panel	41
	6.2		ng parameters	
		6.2.1	Displaying the set temperature for the interior of the device	
		6.2.2	Displaying the offset for the interior temperature	
		6.2.3	Displaying the alarm limits for the interior temperature	
		6.2.4	Displaying the alarm delay time	
		6.2.5	Displaying the ambient temperature	
		6.2.6	Displaying the time for cleaning the air filter	
	6.3		ng with a lock code	
	0.0	6.3.1	Unlocking and locking the device	
		6.3.2	Activating and changing the lock code	
		6.3.3	Deactivating the lock code	
	6.4		mming parameters	
	0.4	6.4.1	Setting the set temperature value for the interior	
		6.4.2	Setting an offset for the set temperature in the interior	
		6.4.3	Setting alarm limits	
		6.4.4	Setting an alarm delay time	
		6.4.5	Setting the interval for cleaning the air filter	
		6.4.6	Confirming the message for cleaning the air filter	
		0.4.0	Commining the message for cleaning the air litter	40
7	Main			
	7.1		e schedule	
	7.2		ting the device	
	7.3	Cleanin	ng and decontamination	50
		7.3.1	Cleaning the device	
		7.3.2	Cleaning and disinfecting the control panel	50
		7.3.3	Cleaning the seals	
		7.3.4	Cleaning the air filter and the air intake grille	51
		7.3.5	Decontaminating the interior and the inner shelves	52
	7.4	Fuses.		53

	7.5	Checkir	ng the alarm	53
		7.5.1	Checking indicator lights and the signal tone	53
		7.5.2	Checking the alarm in the case of a mains/power outage	53
	7.6	Safety o	checklist	53
		•		
8	Troub	leshooti	ing	55
	8.1	General	l errors	55
		8.1.1	Outer door	55
		8.1.2	Air filter	
	8.2	Softwar	e error messages	
	8.3		power outage	
	8.4		up of the interior	
			,	
9	Trans	port, sto	orage and disposal	57
•	9.1	-	missioning	
	9.2		ort	
	· ·-	9.2.1	Preparing the device for transport	
		9.2.2	Transporting the device	
	9.3		nt	
	7.5	9.3.1	Shipping regulations	
		9.3.2	Shipping the device.	
	9.4		il	
	7.4	Dispusa		01
10	Tachr	nical date	a	63
10	10.1		supply	
	10.1		it conditions	
	10.2		Operation	
	10.3		ions	
	10.5	10.3.1	External dimensions	
		10.3.1		
			Internal dimensions.	
	10.4	10.3.3	Packing dimensions	
	10.4	-		
	10.5		evel	
	10.6		res	
	10.7	•	ature control	
			Temperature range	
		10.7.2	Times for cooling and heating the interior	
		10.7.3	Cooling of the refrigeration cycle	
	400	10.7.4	Refrigerant	
	10.8		nal specifications	
		10.8.1	Capacity and carrying capacity	
		10.8.2	Materials	68
11	Order	-	rmation	
	11.1		ories	
		11.1.1	Back-up systems	69
		11.1.2	Chart recorder	
		11.1.3	Racks for devices with 3 compartments	70
		11.1.4	Racks for the compartments 1 – 4 for devices with 5 compartments	70

	11.1.5	Racks for compartment 5 for devices with 5 compartments	71
	11.1.6	Cardboard boxes and box dividers	71
	11.1.7	Eppendorf Storage Boxes	72
Index	· • • • • • • • • • • • • • • • • • • •		73
Certif	icates		75

1 Operating instructions

1.1 Using this manual

- ▶ Read this operating manual completely before using the device for the first time. Observe the instructions for use of the accessories where applicable.
- ▶ This operating manual is part of the product. Please keep it in a place that is easily accessible.
- ▶ Enclose this operating manual when transferring the device to third parties.
- ► The current version of the operating manual for all available languages can be found on our webpage www.eppendorf.com/manuals.

1.2 Danger symbols and danger levels

1.2.1 Danger symbols

The safety instructions in this manual have the following danger symbols and danger levels:

	Risk of tipping over	A	Electric shock
	Highly flammable substances		Explosive substances
*	Low temperatures		Biohazard
	Heavy load		Risk of crushing
<u> </u>	Hazard point	兼	Material damage

1.2.2 Danger levels

DANGER	Will lead to severe injuries or death.
WARNING	May lead to severe injuries or death.
CAUTION	May lead to light to moderate injuries.
NOTICE	May lead to material damage.

1.3 Symbols used

Depiction	Meaning
1.	Actions in the specified order
2.	
•	Actions without a specified order
•	List
Text	Display or software texts
0	Additional information

1.4 Version overview

Version	Date	Change
00	April 2017	Created
01	June 2018	Complete revision
02	June 2018	CryoCube F740h included
03	March 2020	F740 removed New handle and new inner door incorporated
04	August 2020	Warning symbol on the device revised Outer dimensions updated

2 Safety

2.1 Intended use

CryoCube ULT freezers are designed to provide an ultra-low temperature environment for storing scientific research materials. They allow for storage of samples at ultra-low temperatures of -50 °C to -86 °C and at a maximum ambient temperature of 32 °C.

All country-specific safety requirements for operating electrical equipment in laboratories must be observed.

2.2 Warnings for intended use



DANGER! Risk of severe injury from tipping the device over during transport

If the device tips over and falls on someone, that person sustains fatal injuries.

- ▶ Transport the device with a sufficient number of helpers.
- ▶ Observe the transport instructions in the operating manual.



DANGER! Risk of severe injury from climbing onto the device

The device cannot carry the weight of a person. If the device tips over and falls on someone, that person sustains fatal injuries.

The device may become damaged.

- ▶ Do not climb onto the device.
- ▶ Do not pull yourself up on the device or the outer door.



WARNING! Risk of explosion

- ▶ Do not operate the device in areas where work with explosive substances is carried out.
- ▶ Do not store explosive or highly reactive substances in the device.
- ▶ Do not use the device to store substances that may generate an explosive atmosphere.
- ▶ Do not store any aerogenic substances in the device, e.g., dry ice.



WARNING! Lethal voltages inside the device.

If you touch any parts which are under high voltage you may experience an electric shock. Electric shocks cause heart injury and respiratory paralysis.

- Ensure that the housing is closed and undamaged.
- ▶ Do not remove the housing.
- Ensure that no liquids can penetrate the device.

Only authorized service staff may open the device.



WARNING! Electric shock due to damage to the device or unsuitable power cable

If you touch a damaged or unsuitable mains/power cord, you may experience an electric shock. Electric shocks cause injuries to the heart, respiratory paralysis and burns.

▶ If the supplied mains/power cord is defective, replace it with a mains/power cord and a plug of the same type.



WARNING! Damage to health due to infectious liquids and pathogenic germs.

- ▶ When handling infectious liquids and pathogenic germs, observe the national regulations, the biosafety level of your laboratory, and the manufacturers' Safety Data Sheets and application notes.
- ▶ Wear your personal protective equipment.
- ► For comprehensive regulations about handling germs or biological material of risk group II or higher, please refer to the "Laboratory Biosafety Manual" (source: World Health Organization, Laboratory Biosafety Manual, in the currently valid version).



CAUTION! Risk of burns from direct contact with cold surfaces.

The temperature inside the device is low. Direct contact with the interior or samples can cause skin burns.

▶ Wear cold protection gloves when loading and unloading the device.

2.2.1 Devices with flammable refrigerant



WARNING! Risk of fire due to escaping flammable refrigerants (R-170 and R-290)

Refrigerant may leak out if a refrigeration cycle is faulty. The refrigerants R-170 and R-290 are flammable and can form explosive mixtures with the ambient air.

- ▶ Ensure adequate ventilation of the location.
- ▶ Observe the regulations of the owner.
- ▶ Do not allow the device to be maintained or repaired by anyone except authorized service technicians. Components may only be replaced with original spare parts of the same type.

2.3 User profile

The device and accessories may only be operated by trained and skilled personnel.

Before using the device, read the operating manual and the instructions for use of the accessories carefully and familiarize yourself with the device's mode of operation.

2.4 Personal protective equipment

Personal protective equipment protects your life and your health.

- ▶ Always wear the personal protective equipment required for the biosafety level and by the laboratory regulations.
- ▶ Always wear protective clothing, protective gloves, and safety boots.
- ▶ If additional protective equipment is required, this is indicated above the respective instruction.

2.5 Information on product liability

In the following cases, the designated protection of the device may be affected. The liability for any resulting damage or personal injury is then transferred to the owner:

- The device is not used in accordance with the operating manual.
- The device is used outside of its intended use.
- The device is used with accessories or consumables that are not recommended by Eppendorf.
- The device is maintained or repaired by persons not authorized by Eppendorf AG.
- The user makes unauthorized changes to the device.

2.6 Maintenance and repairs

Service technicians authorized by Eppendorf AG are appropriately trained and certified by Eppendorf AG.

- ▶ Do not allow the device to be maintained by anyone except service technicians who are authorized by Eppendorf AG.
 - For more information, please contact your Eppendorf partner or visit www.eppendorf.com.
- ▶ Do not allow the device to be maintained by anyone except service technicians who are accredited according to the national and local laws and safety regulations. Service technicians must hold valid certificates.

Australia, Queensland: the legal regulations state that service technicians require a valid gas work license for working on the refrigeration cycle.

Eppendorf AG uses high-quality components for the device which are manufactured especially for this purpose. These components ensure the safe function of the device. Eppendorf AG provides original spare parts for the service and repair of the device.

▶ Components may only be replaced by original spare parts of the same type.

2.7 Electromagnetic compatibility

2.7.1 Europe

This is a class A product. In a domestic environment this product may cause radio interference in which case the user may be required to take adequate measures.

2.7.2 U.S.A.

Any modification or changes made to this device, unless explicitly approved by Eppendorf, will invalidate the authorization of this device. Operation of an unauthorized device is prohibited under Section 302 of the Communications Act of 1934, as amended, and Subpart I of Part 2 of Chapter 47 of the Code of Federal Regulations.

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the operating manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

2.8 Warning symbols on the device

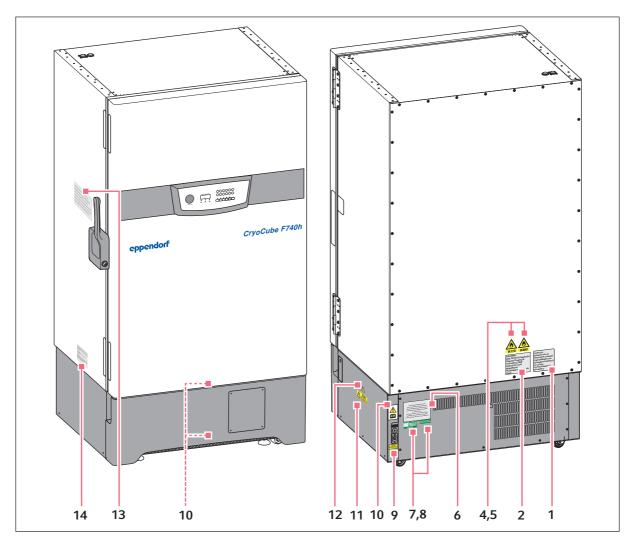


Fig. 2-1: Outside warning symbols

	Warning symbol	Meaning
1	ATTENTION- RISQUE DE FEU OU D'EXPLOSION. ELIMINER CONFORMEMENT AUX REGLEMENTS FEDERAUX OU LOCAUX. LE FRIGORIGENE EST INFLAMABLE	Notice Risk of fire or explosion. Dispose of the device in accordance with federal or local laws and regulations. The device contains flammable refrigerant.
2	CAUTION - Risk Of Fire Or Explosion Dispose Properly In Accordance With Federal Or Local Regulations. Flammable Refrigerants Used.	Notice Risk of fire or explosion. Dispose of the device in accordance with federal or local laws and regulations. The device contains flammable refrigerant.
4	R170	Danger due to flammable refrigerant R-170.
5	R290	Danger due to flammable refrigerant R-290.
6	THIS EPPENDORF FREEZER IS FITTED WITH A CASCADE REFRIGERATION SYSTEM. SERVICE AND REPAIRS MUST BE CARRIED OUT BY A REFRIGERATION SPECIALIST APPROVED BY YOUR SUPPLIER. ANY REPAIRS CARRIED OUT BY UNAUTHORISED ENGINEERS COULD CAUSE SERIOUS DAMAGE TO THE SYSTEM AND MAY AFFECT YOUR WARRANTY.	The device may only be serviced and repaired by a qualified refrigeration specialist who has been authorized by Eppendorf AG. If the device is serviced or repaired by an unauthorized person, liability on the part of Eppendorf AG shall cease immediately.

	Warning symbol	Meaning
7	DATE/INITIALS APPL/REF NO. NEXT TEST DATE PASSED	The device has passed the electrical safety test.
8	RoHS Compliant Directive 2011/65/EU	The device complies with the Restriction of Hazardous Substances (RoHS) Directive 2011/65/EU.
9	WARNING THIS EQUIPMENT MUST BE EARTHED	Only connect the device to a mains/power connection with PE conductor.
10		Notice of a hazard point. Read the operating manual.
11	Danger Disconnect the mains supply before removing this cover	Risk of electric shock. Disconnect the mains/power cord from the voltage supply before removing the cover.
12	A	Risk of electric shock.

Warning symbol Meaning 13 Notice of a hazard point. Read the operating manual: ICE FORMATION ON SEAL! ICE FORMATION ON SEAL! After prolonged or frequent opening of the freezer, humidity or ice may form on the door seal and the case frame, Ice formation may impair the function of the high-efficiency door seal, • Wipe off humidity before closing the door, • Remove Ice from the sealing surfaces. Risk of crushing injuries when closing the outer door. Ice formation on the seal VACUUM INSIDE FREEZER! After closing the outer door, a vacuum can occur inside the freezer. The outer door can no longer be opened. To speed up pressure compensation, press the *auto* verif button. Wait until pressure compensation has taken place, then operate the handle. Opening the device often or leaving it open for a longer period of time will lead to moisture or ice forming on the door seal and on the housing. Ice can impair the function of the door seal. • Wipe off any moisture before closing the door. • Remove any ice from the surface of the seal. Negative pressure in the interior A negative pressure can be generated after closing the outer door. The outer door can no longer be opened. • To speed up pressure equalization, press the *auto vent* • Wait for the pressure to equalize. Use the door handle only after this has occurred. 14 The device complies with FCC Rules Part 15. THIS DEVICE COMPLIES WITH PART15 OF THE FCC RULES. OPERATION IS SUBJECT TO THE FOLLOWING CONDITIONS: The device may be operated under the following (1) THIS DEVICE MAY NOT CAUSE HARMFUL INTERFERENCE conditions: • This device may not cause harmful interference. (2) THIS DEVICE MUST ACCEPT ANY INTERFERENCE RECEIVED, INCLUDING INTERFERENCE THAT MAY CAUSE UNDESIRED OPERATION • This device must accept any interference received. This includes interference that may cause undesired operation.

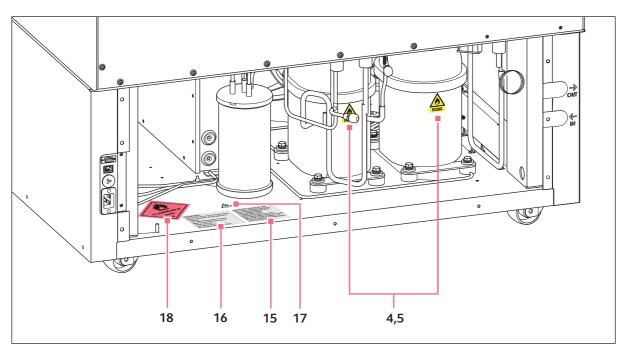


Fig. 2-2: Warning symbols on the rear of the opened device

The warning symbols inside the device can only be accessed by authorized service technicians.

	Warning symbol	Meaning
15	DANGER-RISQUE DE FEU OU D'EXPLOSION. LE FRIFORIGENE EST INFLAMMABLE. CONFIER LES REPARATIONS A UN TECHNICIEN SPECIALISE. NE PAS PERFORER LA TUBULURE CONTENANT LE FRIGORIGENE.	Danger Risk of fire or explosion The device contains flammable refrigerant. Only qualified service technicians are allowed to repair the device. Do not puncture the refrigerant tubing.
16	DANGER - Risk Of Fire Or Explosion Flammable Refrigerant Used. To Be Repaired Only By Trained Service Personnel. Do Not Puncture Refrigerant Tubing.	Danger Risk of fire or explosion The device contains flammable refrigerant. Only qualified service technicians are allowed to repair the device. Do not puncture the refrigerant tubing.

	Warning symbol	Meaning
17		Danger due to flammable materials
18	FLAMMABLE GAS 2	Danger due to flammable gas, class 2

3 Product description

3.1 Product overview

3.1.1 Front view

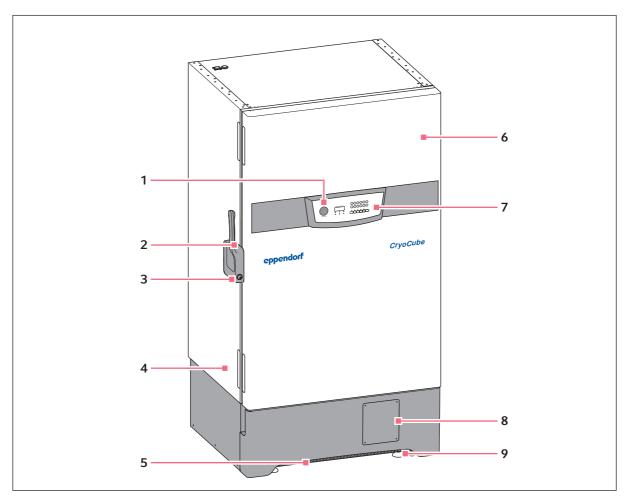


Fig. 3-1: Front view of a model with the door handle mounted on the left side

- 1 *auto vent* valve
 Automatic pressure compensation
- 2 Door handle
- 3 Mechanical lock
- 4 Name plate
- 5 Air filter

- 6 Outer door
- 7 Control panel
 Display and operating controls
- 8 Space for an optional chart recorder
- 9 Leveling feet

3.1.2 Side view

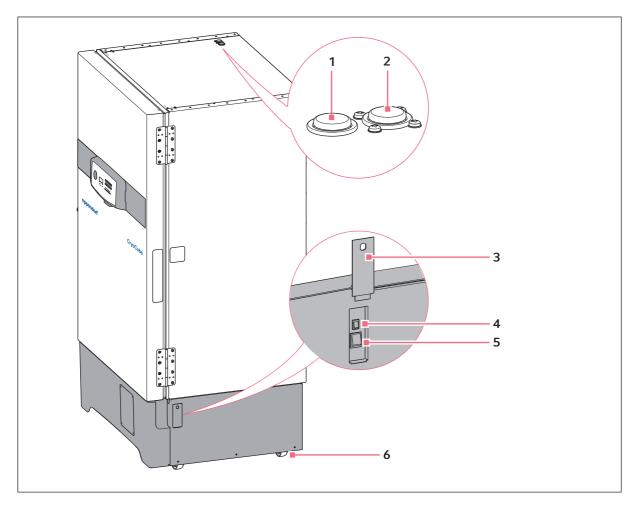


Fig. 3-2: Side view

1 Access port

For external sensors

2 Access port

For external sensors or an optional back-up system

3 Switch locking plate

Lockable plate cover for the mains/power switch and the battery switch

4 Battery switch

For activating the back-up circuit

- 5 Mains/power switch
- 6 Heavy-duty castors

3.1.3 Internal view

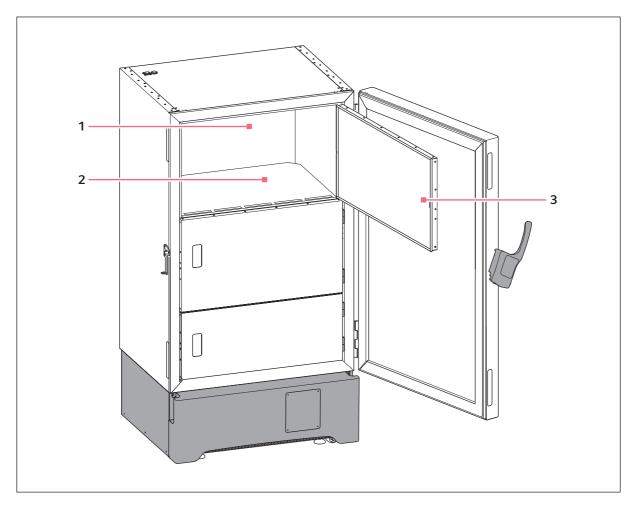


Fig. 3-3: Internal view of a model with three inner doors

1 Inner compartment

3 Inner door with snap latch

2 Shelf

3.1.4 Interfaces

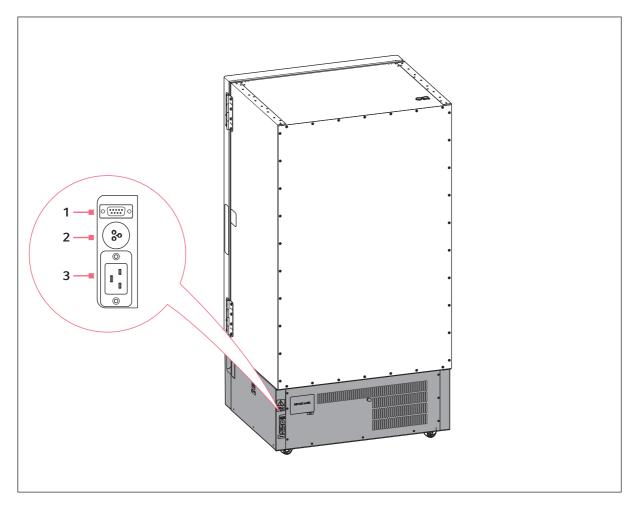


Fig. 3-4: Rear view

- 1 Serial interface
 Connection for internal use
- 2 Remote alarm interface
 Connection to a building management system
- 3 Mains/power connection Connection for the mains/power cord

Only connect devices that comply with the IEC 60950 (UL 60590) standards to the interfaces.

3.2 Features

The CryoCube is a ULT freezer for storing biological samples.

The device has a two-stage cascade refrigeration system with two closed refrigeration cycles. The refrigeration cycles are air-cooled.

The device is operated using the control panel. The control panel consists of a display, softkeys and indicator lights.

The device communicates with external systems over the interfaces. Alarms and other parameters are transmitted.

The outer door is locked and unlocked manually.

The device features automatic pressure compensation. Automatic pressure compensation occurs as soon as the outer door is closed. Afterwards, the outer door can be opened again. Pressure compensation can also be triggered with the *auto vent* valve.

Each inner door is equipped with a seal. With the inner doors closed, cold air loss is minimized and ice formation inside the device is reduced.

The devices are equipped with 2 or 4 shelves. The number and position of the shelves cannot be changed.

The device is equipped with a switch locking plate. The mains/power switch and the battery switch are located behind this lockable cover.

An air filter is located below the outer door. The air filter protects the condenser and the assemblies behind it against contamination and dirt.

The heavy-duty castors serve to transport the device safely to its place of installation. Leveling feet ensure the stability and horizontal alignment of the device.

3.3 Models

Name	Door handle	Number of compartments	Cooling
CryoCube F740h	Left	3	Air cooling
CryoCube F740h	Left	5	Air cooling
CryoCube F740h	Right	3	Air cooling
CryoCube F740h	Right	5	Air cooling

3.4 Alarms

When a safety-relevant situation has been reached, the device triggers an alarm. The user must immediately eliminate the cause of the alarm.

Alarms consist of a signal tone on site, an indicator light on the control panel, an alarm message on the control panel and forwarding of the alarm to a building management system. As soon as the cause of the alarm has been eliminated, all alarm signals will stop.

Hazard messages can be configured (see *Setting alarm limits on p. 46*), (see *Setting an alarm delay time on p. 46*).

Alarm: Interior temperature

- The temperature in the interior exceeds the alarm limit for the maximum or minimum temperature in the interior.
- The alarm is triggered when the delay time has elapsed. The delay time can be defined.
- The signal tone sounds on the device.
- The **temp alarm** indicator light lights up.
- The alarm is forwarded to a building management system.
- The alarm cannot be deactivated.

Alarm: Mains/power outage

- The mains/power supply to the device is interrupted. The battery back-up circuit is switched on and triggers the alarm.
- The signal tone sounds on the device.
- The **power fail** indicator light lights up.
- The display shows the interior temperature and flashes at intervals of 10 s.
- The alarm is forwarded to a building management system.
- The alarm cannot be deactivated.

Alarm: Battery

- The battery voltage is too low.
- The **battery-low** indicator light lights up.
- The alarm cannot be deactivated.

Alarm: System error

- The signal tone sounds on the device.
- The fault indicator light lights up.
- The alarm cannot be deactivated.

Alarm: Clean the air filter

- The servicing date for the air filter has been reached. The air filter has to be cleaned.
- The filter-clean indicator light flashes.
- The alarm cannot be deactivated.

3.5 Delivery package

3.5.1 Device and accessories

Quantity	Description
1	ULT upright freezer
1 or 2	Mains/power cord (quantity depends on country)
1	Safety clamp for mains/power cord
2	Key for switch locking plate
2	Key for outer door
2	Anti-slip pads
1	Plug for connection to the building management system
1	Allen key

3.5.2 Documents

Quantity	Description
1	Operating manual
1	Unpacking instructions
1	Certificate of conformity

3.6 Accessories

Optional accessories can be ordered separately. Information regarding accessories is available on our website: www.eppendorf.com.

3.6.1 Back-up systems

In the event of a mains/power outage, the battery-powered back-up system will start and cool the interior for a limited period of time. The back-up system is connected through the access port.

The following back-up systems are available:

- $\rm CO_2$ back-up system for temperatures from -60 °C to -70 °C.
- LN₂ back-up system for temperatures down to -85 °C.

3.6.2 Chart recorder

The chart recorder records the interior temperature on a disk over a period of 7 days. The port for connection of the chart recorder is available in the device.

Pens and disks for chart recorders are available.

3.6.3 Racks for ULT upright freezers

Racks serve to store and sort samples in boxes, microplates and deepwell plates.

Racks are placed on the inner shelves inside the device. The racks are used to store boxes. Stainless steel racks provide space for $136 \text{ mm} \times 136 \text{ mm}$ boxes. Aluminum racks can be used to store boxes up to $133 \text{ mm} \times 133 \text{ mm}$.

Racks are available with drawers or with open sides. The design of the racks ensures an even temperature in the entire rack.

3.6.4 Cardboard boxes and box dividers

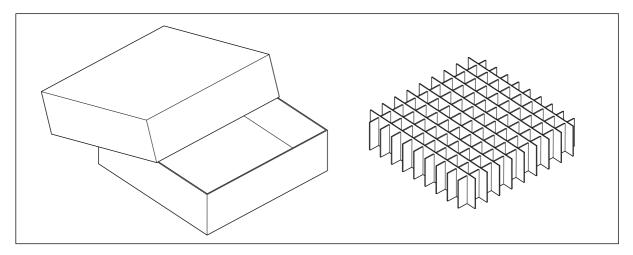


Fig. 3-5: Cardboard box and box divider

Cardboard boxes serve to store samples in tubes at temperatures down to -86 °C. Cardboard boxes have a waterproof coating.

To sort your samples, you can insert box dividers into the cardboard boxes. Eppendorf AG cardboard boxes and box dividers are compatible with each other.

3.6.5 Eppendorf Storage Box

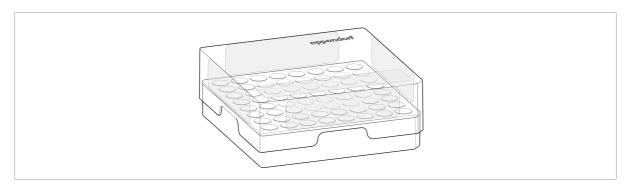


Fig. 3-6: Eppendorf Storage Box

Eppendorf Storage Boxes serve to store samples in tubes at temperatures down to -86 °C.

Eppendorf Storage Boxes are made of polypropylene (PP) and are autoclavable.

Product description CryoCube® F740h English (EN)

28

4 Installation

4.1 Selecting the location

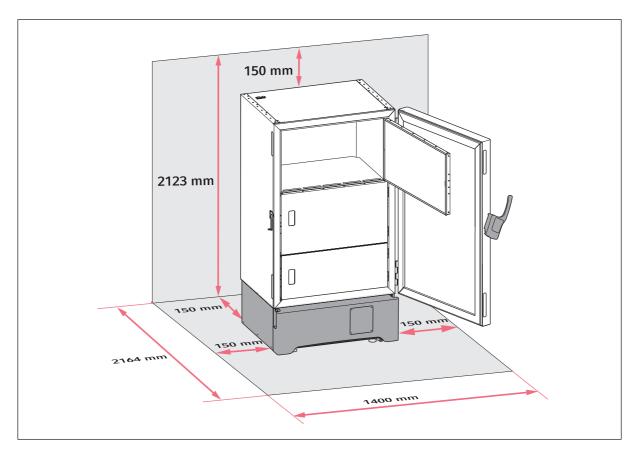


Fig. 4-1: Footprint

Information on ambient conditions, dimensions and weights can be found in the technical data.

Location in general

- The ambient conditions match the specifications set out in the Technical data chapter.
- The location is well ventilated or air-conditioned.
- The location is not next to heat sources.
- The location is protected against sparks and open fire.
- The floor space corresponds to the technical data.
- The floor is level, vibration-free and designed for the weight of the device.

Electrical connection

- Mains/power connection in accordance with the name plate
- The mains/power switch of the device and the disconnecting device of the power system circuit (e.g., residual current circuit breaker) are accessible during operation.



Contact your safety officer for information on further requirements when installing the device.

4.2 Preparing installation

4.2.1 Unpacking the device

- 1. Check the packaging for damage.
- 2. Unpack the device in accordance with the unpacking instructions.

4.2.2 Checking the delivery

- 1. Check the delivery for completeness.
- 2. Check the device and accessories for transport damage.
- 3. Do not commission the device if the packing or the device is damaged. Contact Eppendorf AG customer service or your Eppendorf partner.

4.2.3 Transporting the device to the location

Personal protective equipment

• Protective clothing, protective gloves, safety shoes

Prerequisites

- The location meets the requirements.
- ▶ Transport the device to the location (see *Transport on p. 57*).

4.2.4 Setting up the device



CAUTION! Risk of cut injuries due to sharp-edged components

When aligning the leveling feet, cut injuries can result from sharp-edged components on the bottom of the device.


▶ Wear protective gloves when setting up the device.

Personal protective equipment


• Protective clothing, protective gloves, safety shoes

Prerequisites

• The device is in its intended position.

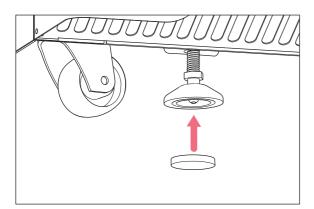

1. Attach anti-slipping pads to the underside of the leveling feet.

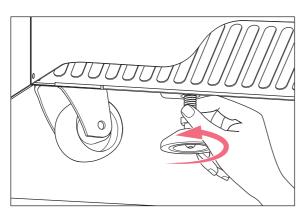

2. Rotate the leveling feet down.

3. Remove the adhesive tape from the air intake grille.

4.3 Connecting the device to the voltage supply



WARNING! Risk due to incorrect voltage supply.

- ▶ Only connect the device to voltage sources which correspond with the electrical requirements on the name plate.
- ▶ Only use earth/grounded sockets with a protective earth (PE) conductor.
- ▶ Only use the mains/power cord supplied.


Prerequisites

- Mains/power connection in accordance with the name plate
- 1. If several mains/power cords are included, select the mains/power cord according to the mains/power supply voltage.
- 2. Connect the mains/power cord at the rear of the device.
- 3. Fasten the safety clamp.

If the device is moved, the mains/power cord cannot be pulled out of the device.

4.4 Connecting the device to external systems

4.4.1 Remote alarm interface

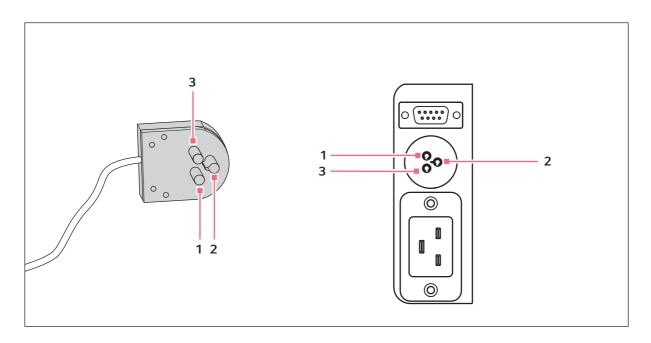

- 1 Pin 1 and socket 1
- 2 Pin 2 and socket 2

3 Pin 3 and socket 3

You can connect the device to a building management system via the remote alarm interface.

The following alarms are forwarded to the building management system:

- Alarm in the case of a mains/power outage
- Alarm that the temperature inside the device is too high
- Alarm that the temperature inside the device is too low

The plug is included in the delivery package. Connections must have double or reinforced insulation as described in DIN EN 61010-1.

4.4.2 RS-485 interface

The RS-485 interface is reserved for internal use.

4.5 Switching the device on



WARNING! Risk of electric shock due to damage to the device or to the mains/power cord.

- ▶ Only switch on the device if the device and the mains/power cord are undamaged.
- ▶ Only operate devices which have been installed or repaired properly.
- ▶ In the event of danger, disconnect the device from the mains/power supply voltage.

 Disconnect the mains/power plug from the device or the earth/grounded socket. Use the isolating device intended for this purpose (e.g., the emergency switch in the laboratory).



NOTICE! Damage to electronic components due to condensation.

Condensate may form in the device when it has been transported from a cool environment to a warmer environment.

▶ After installing the device, wait for at least 6 h. Only then connect the device to the mains/ power line.



NOTICE! Improper door sealing due to ice

Humidity inside the device causes ice formation. Ice causes damage to the seals of the inner and outer doors.

- 1. Dry the device completely, especially all seals.
- 2. Switch on the device afterwards.

4.5.1 Enabling the back-up circuit

The back-up circuit is battery-powered. In the event of a mains/power outage, the battery will supply power to the control panel and the alarm for 72 hours.

Tools and auxiliary equipment

• Key for the switch locking plate

Prerequisites

- The device has been installed and connected according to the operating manual.
- The device has been acclimatized for at least 6 h.
- 1. Unlock the switch locking plate and remove the cover.
- 2. Switch on the battery switch.

The back-up circuit is activated.

- An alarm is triggered in the event of a mains/power outage.
- In the event of a mains/power outage, power will still be supplied to the control panel.
- The software settings are saved in the event of a mains/power outage.
- The battery is charged with mains power. The battery is fully charged after approx. 24 hours.
- 3. Fit the cover and lock it.

4.5.2 Switching the device on at the mains/power switch

Tools and auxiliary equipment

• Key for the switch locking plate

Prerequisites

- The device has been installed and connected according to the operating manual.
- The device has been acclimatized for at least 6 h.
- The interior, seals, doors and lids are dry.
- 1. Unlock the switch locking plate and remove the cover.
- 2. Switch on the mains/power switch.
 - The display shows the software version number.
 - The compressor starts running after a short time delay.
- 3. Fit the cover and lock it.

4.6 Basic device settings

To commission the device, set the following values. Further settings are described in the Software chapter.

- 1. Set the desired temperature for the interior of the device.
- 2. Set the temperature offset.
- 3. Set alarm limits.
- 4. Set a delay time for the alarm.
- 5. Check the alarms.

Installation CryoCube® F740h English (EN)

36

5 Operation

5.1 Opening the outer door




CAUTION! Hand injuries

There are moving parts on the inside of the door handle.

▶ Do not touch the inside of the door handle.

Prerequisites

• Pressure compensation has finished.

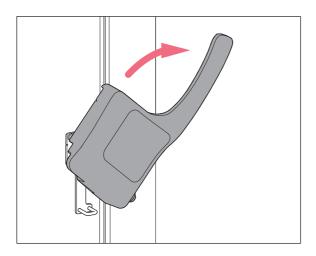

- 1. Unlock and remove the padlock, where applicable.
- 2. Pull the door handle forward and down until it stops.
- 3. To open the outer door, pull the door handle forward.



If the outer door will not open, press the *auto vent* button to accelerate pressure equalization.

5.2 Loading the device



CAUTION! Risk of head injury due to open inner door

If the upper inner doors are opened you can hit your head on the inner doors.

- ▶ Only open one inner door at a time.
- ▶ Immediately close the inner door after completing your work.



CAUTION! Risk of fingers being crushed when closing the inner doors.

Incorrect handling of the inner doors can cause crushing injuries.

- ▶ Only grasp the inner doors by the handle.
- ▶ Always open only one inner door.



NOTICE! Longer pull-down time because the device is loaded too early

The pull-down time is the time needed for the device to cool the interior from the ambient temperature to the set temperature.

If you load the device during the cooling phase, the pull-down time will be longer. The pull-down time specified in the technical data cannot be achieved.

- 1. Allow the device to cool down from ambient temperature to the set temperature.
- 2. Place the samples in the device after the device has reached the set temperature.



The interior temperature of the device increases when loading it:

- · Outer and inner doors are open.
- The sample temperature differs from the interior temperature.
- ▶ To minimize the temperature increase in the interior, load the device step by step.

Prerequisites

- Racks and accessories have been placed in the compartments.
- Device, racks and accessories have reached the set temperature.
- 1. Open the outer door.
- 2. Open the inner door of the compartment in which you want to place the samples.
- 3. Place the samples in the device.

 Information on the maximum carrying capacity of the inner shelves can be found in the technical data.
- 4. Close the inner door.
- 5. Close the outer door.

5.3 Locking the outer door



CAUTION! Risk of hand being crushed when closing the outer door

- ▶ Do not place your fingers between the device and the outer door.
- ▶ Lock the door handle slowly and carefully.




NOTICE! Damage to the door handle due to incorrect closing of the outer door.

Closing the outer door while the door handle is in the upright position damages the door handle.

- 1. Pull the door handle forward and down first.
- 2. Then close the outer door.



If the door handle is not properly engaged and there is negative pressure in the interior, the outer door is closed. However, as soon as the negative pressure is compensated, the outer door opens again.

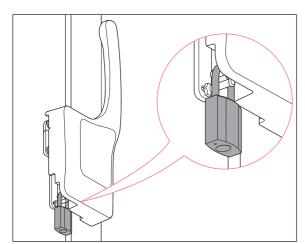

- 1. Press the door handle forward and down.
- 2. Close the outer door.
- 3. Lock the outer door. To do so, press the door handle up.
 - Automatic pressure compensation takes place as soon as the outer door is closed.
- 4. Check if the door handle is engaged in locked position.
- 5. Mount the padlock and lock it, where applicable.

5.4 Pressure compensation

If you leave the outer door open for a while, the temperature in the interior will increase. After closing the outer door, the air in the interior will cool down and the atmospheric pressure will decrease. Negative pressure may occur in the device. If negative pressure occurs, the outer door can no longer be opened.

Pressure compensation takes place automatically to reestablish ambient pressure in the device. Pressure compensation starts as soon as the outer door is closed.

➤ To speed up pressure compensation, press the *auto vent* valve.

With the *auto vent* valve, pressure compensation takes 1 min – 2 min.

5.5 Switching off the device



WARNING! Risk of electric shock due to damage to the device or to the mains/power cord.

- ▶ Only switch on the device if the device and the mains/power cord are undamaged.
- ▶ Only operate devices which have been installed or repaired properly.
- ▶ In the event of danger, disconnect the device from the mains/power supply voltage.

 Disconnect the mains/power plug from the device or the earth/grounded socket. Use the isolating device intended for this purpose (e.g., the emergency switch in the laboratory).

5.5.1 Disabling the back-up circuit

Tools and auxiliary equipment

- · Key for the switch locking plate
- 1. Unlock the switch locking plate and remove the cover.
- 2. Switch off the battery switch.

The back-up circuit is disabled.

- No alarm is triggered in the event of a mains/power outage.
- During a mains/power outage, there will be no power supply to the control panel.
- · The battery is not charged.

5.5.2 Disconnecting the device from the voltage supply

Tools and auxiliary equipment

- Key for the switch locking plate
- 1. Unlock the switch locking plate and remove the cover.
- 2. Switch off the battery switch.

The back-up circuit is disabled.

- No alarm is triggered in the event of a mains/power outage.
- During a mains/power outage, there will be no power supply to the control panel.
- The battery is not charged.
- 3. Switch off the mains/power switch.

6 Software

6.1 Overview of the control panel

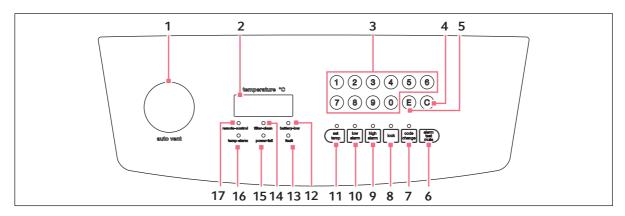


Fig. 6-1: Control panel

1 auto vent valve

Automatic pressure compensation

2 Display

During operation the display shows the actual temperature in the interior.

3 Numerical keys

Access functions. Enter values.

4 Ckey

Delete entry.

5 E key

Confirm your entry.

6 alarm test mute softkey

Disable the signal tone. Test the signal tone.

7 code change softkey

Change the lock code.

8 lock softkey

Lock or unlock control panel.

9 high alarm softkey

Display and adjust the alarm limit for the maximum interior temperature.

10 low alarm softkey

Display and adjust the alarm limit for the minimum interior temperature.

11 set temp softkey

Display and adjust the set interior temperature.

12 battery-low indicator light

Lights up or flashes if the voltage in the back-up circuit is too low.

Lights up if no mains/power supply voltage is available.

13 fault indicator light

Lights up if a system error has occurred.

14 filter clean indicator light

Lights up if the air filter needs to be cleaned.

15 power fail indicator light

Flashes if no mains/power supply voltage is supplied to the device.

The signal tone sounds while the indicator light is flashing.

16 temp-alarm indicator light

Lights up if an alarm limit for the interior temperature has been exceeded.

17 remote control indicator light

Lights up if the device is controlled by means of a computer.

6.2 Checking parameters

6.2.1 Displaying the set temperature for the interior of the device

▶ Press the **set temp** softkey.

The set temperature for the interior of the device is displayed. The factory setting is -80 °C.

6.2.2 Displaying the offset for the interior temperature

▶ Press the C softkey.

The offset for the interior temperature is displayed.

6.2.3 Displaying the alarm limits for the interior temperature

- ▶ To display the alarm limit for the maximum interior temperature, press the **high alarm** softkey.
- ▶ To display the alarm limit for the minimum interior temperature, press the **low alarm** softkey. The alarm limit is displayed.

6.2.4 Displaying the alarm delay time

- ▶ To display the delay time for the on-site alarm, press the **8** softkey.
- ► To display the delay time for the remote alarm, press the **9** softkey. The delay time is displayed.

6.2.5 Displaying the ambient temperature

▶ Press the **0** softkey.

The ambient temperature is displayed.

6.2.6 Displaying the time for cleaning the air filter

Press the 7 softkey.

The display shows the time remaining until the next cleaning is due in months.

6.3 Working with a lock code

To prevent unauthorized programming of the device, you can use a four-digit lock code.

6.3.1 Unlocking and locking the device

Unlocking the device

Prerequisites

- The lock code is activated.
- 1. Press the **lock** softkey.

If the lock indicator light flashes, a lock code has been set.

2. Enter the lock code.

The actual temperature is displayed.

The lock indicator light glows. The device is in programming mode. The parameters can be changed.

Locking the device

Prerequisites

- The lock code is activated.
- 3. After programming is completed, press the **lock** softkey.

The **lock** indicator light goes out. The device is no longer in programming mode. The parameters are saved.



If you press a softkey, e.g., **set temp**, while the **lock** indicator light is flashing, ---- appears on the display. The control panel is locked.

6.3.2 Activating and changing the lock code



NOTICE! No software access due to lost lock code

If you lose the lock code, reprogramming of the device will no longer be possible. The lock code will have to be reset by an authorized service technician.

▶ Keep the lock code in a safe place.

In the factory setting, the lock code is deactivated. To activate or change the lock code, proceed as follows.

- 1. Press the **lock** softkey.
- 2. When the lock indicator light flashes, enter the lock code.

The lock indicator light glows. The device is in programming mode. The parameters can be changed.

3. Press the **code change** softkey.

The **code change** indicator light flashes. The display is empty.

4. Enter a lock code using the numerical keys.

The lock code is displayed.

- 5. Check the lock code on the display.
- 6. To delete the entry, press softkey C.

7. Confirm the entry. To do so, press softkey **E**.

The code change indicator light goes out.

The new lock code is active.

8. Exit programming mode. To do so, press the lock softkey.

The **lock** indicator light goes out. The device is no longer in programming mode. The parameters are saved.

6.3.3 Deactivating the lock code

To deactivate the lock code, set it to **0000**.

1. Press the lock softkey.

The lock indicator light flashes.

2. Enter the current lock code.

The actual temperature is displayed.

The lock indicator light glows. The device is in programming mode. The parameters can be changed.

3. Press the code change softkey.

The code change indicator light flashes. The display is empty.

4. Enter the lock code **0000** using the numerical keys.

Entering the lock code **0000** will deactivate the lock code.

The lock code 0000 is displayed.

- 5. Check the lock code on the display.
- 6. To delete the entry, press softkey C.
- 7. Confirm the entry. To do so, press softkey **E**.

The **code change** indicator light goes out.

The new lock code is deactivated.

8. Exit programming mode. To do so, press the lock softkey.

The **lock** indicator light goes out. The device is no longer in programming mode. The parameters are saved.

6.4 Programming parameters

6.4.1 Setting the set temperature value for the interior

You can set the set temperature for the interior of the device to a range from -50 °C to -86 °C.

Prerequisites

- The device is not protected by a lock code.
- 1. Press the lock softkey.

The lock indicator light lights up. The device is in programming mode. The parameters can be changed.

2. Press the **set temp** softkey.

The **set temp** indicator light flashes. *0* is displayed.

3. Enter the set temperature with the numerical keys.

The set temperature is displayed. The set temperature is automatically displayed as a negative value.

- 4. To delete the entry, press the **C** softkey.
- 5. Confirm the entry. To do so, press the **E** softkey.

The **set temp** indicator light goes out.

6. Exit programming mode. To do so, press the lock softkey.

The lock indicator light goes out. The parameters are saved.

6.4.2 Setting an offset for the set temperature in the interior

You can set an offset of 0 °C to -5 °C for the set temperature value in the interior. The offset is added to the set temperature. The device temperature cannot fall below -86 °C.

Prerequisites

- The device is not protected by a lock code.
- 1. Press the **lock** softkey.

The lock indicator light lights up. The device is in programming mode. The parameters can be changed.

- 2. Press the C softkey.
- 3. Enter the offset using the numerical keys.

The offset is displayed.

To deactivate the offset, enter the value 0.

- 4. To delete the entry, press the **C** softkey.
- 5. Confirm the entry. To do so, press the **E** softkey.
- 6. Exit programming mode. To do so, press the **lock** softkey.

The **lock** indicator light goes out. The parameters are saved.

6.4.3 Setting alarm limits

You can set alarm limits for the interior temperature. If the interior temperature exceeds an alarm limit, an alarm will be triggered.

	Minimum value	Maximum value
Alarm limit for the minimum temperature in the interior	-91 °C	5 °C below the set temperature
Alarm limit for the maximum temperature in the interior	5 °C above the set temperature	-10 °C

Prerequisites

- The device is not protected by a lock code.
- 1. Press the lock softkey.

The lock indicator light lights up. The device is in programming mode. The parameters can be changed.

- 2. To set the alarm limit for the maximum interior temperature, press the **high alarm** softkey.
 - The high alarm indicator light flashes. The display shows $\theta.$
- 3. To set the alarm limit for the minimum interior temperature, press the **low alarm** softkey. The **low alarm** indicator light flashes. *0* is displayed.
- 4. Enter the alarm limit using the numerical keys.

The display shows the alarm limit.

- 5. To delete the entry, press the **C** softkey.
- 6. Confirm the entry. To do so, press the **E** softkey.

The high alarm indicator light goes out.

7. Exit programming mode. To do so, press the **lock** softkey.

The lock indicator light goes out. The parameters are saved.

6.4.4 Setting an alarm delay time

You can set a delay time for the "Interior temperature too high" and "Interior temperature too low" alarms. The delay time can be set for the on-site alarm and the remote alarm.

	Minimum value	Maximum value	Factory setting
On-site alarm	0 min	40 min	30 min
Remote alarm	0 min	40 min	30 min

If you set the delay time to 0 min, the delay time will automatically be set to 15 s.

Prerequisites

- The device is not protected by a lock code.
- 1. Press the **lock** softkey.

The lock indicator light lights up. The device is in programming mode. The parameters can be changed.

- 2. To set the delay for the on-site alarm, press the 8 softkey.
- 3. To set the delay for the remote alarm, press the **9** softkey.

The PP value is displayed.

4. Enter the delay time using the numerical keys.

The delay time is displayed.

- 5. To delete the entry, press the **C** softkey.
- 6. Confirm the entry. To do so, press the **E** softkey.
 - --- is displayed. The value is saved.
- 7. If the entered value is outside the thresholds, the message *EE* is displayed. Repeat the entry.
- 8. Exit programming mode. To do so, press the lock softkey.

The lock indicator light goes out. The parameters are saved.

6.4.5 Setting the interval for cleaning the air filter

You can set the interval for cleaning the air filter to a range from 0 months to 12 months. The factory setting is 3 months.

Prerequisites

- The device is not protected by a lock code.
- 1. Press the lock softkey.

The lock indicator light lights up. The device is in programming mode. The parameters can be changed.

- 2. Press the alarm test/mute softkey.
- 3. Enter the interval for the air filter check.

The display shows the interval.

To deactivate the function, enter the value 0.

- 4. To delete the entry, press the **C** softkey.
- 5. Confirm the entry. To do so, press the **E** softkey.
- 6. Exit programming mode. To do so, press the lock softkey.

The lock indicator light goes out. The parameters are saved.

6.4.6 Confirming the message for cleaning the air filter

When the interval for cleaning the air filter has elapsed, the **filter clean** indicator light will start to flash.

- 1. Clean the air filter (see Cleaning the air filter and the air intake grille on p. 51).
- 2. Press the **7** softkey for 10 seconds.

The filter clean indicator light goes out.

The display shows the time remaining until the next cleaning is due in months. Time counting restarts from the beginning.



You can also restart the time counter from the beginning without cleaning the air filter.

7 Maintenance

7.1 Service schedule

Service	Service cycle
Defrost the device.	As required
Clean the interior and exterior of the device.	As required
Clean the seals.	Once a month
Clean the air filter and the air intake grille.	Every 3 months under normal ambient conditions. Clean more frequently if the surroundings are very dusty or dirty.

7.2 Defrosting the device



CAUTION! Risk of slipping due to melt water

Puddles may form on the laboratory floor when defrosting the device.

▶ Wipe up melt water immediately.



NOTICE! Risk of device damage due to scraping off ice

Removing ice with a sharp object may damage the device.

▶ Wait until the ice has thawed by itself.



NOTICE! Improper door sealing due to ice

Humidity inside the device causes ice formation. Ice causes damage to the seals of the inner and outer doors.

- 1. Dry the device completely, especially all seals.
- 2. Switch on the device afterwards.

Tools and auxiliary equipment

- Personal protective equipment: Cold protection gloves, protective goggles, dust protection mask
- Material for absorbing the melt water
- "Defrosting device" notice sign

Prerequisites

- The samples have been transferred to another ULT freezer.
- The device is switched off and disconnected from the mains/power line (see p. 40).
- 1. Put up the notice sign.
- 2. Open the outer and inner doors of ULT upright freezers. Open the outer and inner lids of ULT chest freezers.
- 3. Wait until the ice has thawed.
- 4. Wipe up the melt water.
- 5. Dry the device completely, especially all seals.

7.3 Cleaning and decontamination



DANGER! Electric shock.

▶ Switch off the device and disconnect the mains/power plug before commencing any service or cleaning procedures.



NOTICE! Damage from the use of aggressive chemicals.

- ▶ Do not use any aggressive chemicals on the device or its accessories, such as strong and weak bases, strong acids, acetone, formaldehyde, halogenated hydrocarbons or phenol.
- ▶ If the device has been contaminated by aggressive chemicals, clean it immediately using a mild cleaning agent.

7.3.1 Cleaning the device

Tools and auxiliary equipment

- Water
- · Mild cleaning agent
- · Soft, lint-free cloth

Prerequisites

- For cleaning the interior: The device is switched off and disconnected from the mains/power line.
- The device is defrosted.
- 1. If required, lift the inner shelf upwards out of the device.
- 2. Moisten the lint-free cloth with water and cleaning agent.
- 3. Clean the surfaces.

7.3.2 Cleaning and disinfecting the control panel

Tools and auxiliary equipment

- · Laboratory cleaner
- Lint-free cloth
- Disinfectant: Ethanol 70%, sodium hypochlorite solution 1%, Dismozon pur, Hexaquart S, Biozid ZF or another suitable disinfectant
- 1. To lock the control panel, press the **lock** softkey.
- 2. Moisten the lint-free cloth with laboratory cleaner or disinfectant.
- 3. Wipe the control panel with the cloth.
- 4. Unlock the locked control panel.

7.3.3 Cleaning the seals

Tools and auxiliary equipment

- · Dry soft lint-free cloth
- 1. Wipe the seal with a soft, lint-free cloth.
- 2. Wipe the contact surface for the seal with a soft, lint-free cloth.

7.3.4 Cleaning the air filter and the air intake grille



CAUTION! Risk of cut injuries due to sharp-edged components

When working on the air intake grille, cut injuries can be caused by sharp-edged components.

▶ Wear protective gloves to avoid cut injuries.



NOTICE! Refrigeration failure due to blocked air filter

If the air filter is blocked, the refrigerant will not be liquefied. This will damage the compressor.

▶ Regularly check that the air flow into the device is not obstructed.

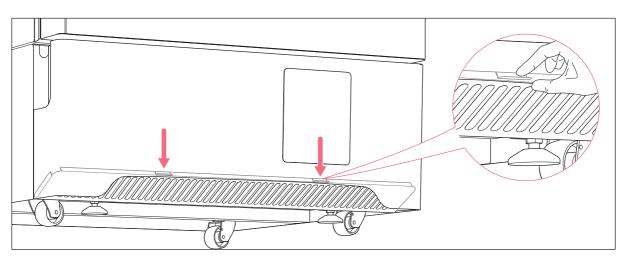


Fig. 7-1: Folding down the air intake grille

An airflow from the surroundings flows through the air filter to the condenser. The airflow carries the heat away from the condenser.

The air filter protects the condenser and the assemblies behind it against contamination and dirt. If the air filter is dirty, not enough air will reach the condenser. The condenser will overheat and the refrigeration cycle may fail.

Personal protective equipment

· Protective gloves

Tools and auxiliary equipment

- · Vacuum cleaner
- · Warm water
- 1. Set the interval for cleaning the air filter in the software.

When the interval has elapsed, the filter-clean indicator light will start to flash.

- 2. Place your fingers into the recesses on the air intake grille. Press the air intake grille downwards. The air intake grille folds down.
- 3. Remove the air filter.
- 4. Clean the air intake grille with the vacuum cleaner.
- 5. Alternatively, clean the air intake grille with a soft brush.
- 6. Remove the coarse dirt from the air filter by vacuuming or tapping it.
- 7. Clean the air filter with warm water.
- 8. Let the air filter dry.
- 9. Insert the air filter.
- 10. Press the air intake grille upwards and close it.

7.3.5 Decontaminating the interior and the inner shelves

The interior and the shelves are made of stainless steel.

Tools and auxiliary equipment

- Decontamination agent consisting of 70 % isopropyl alcohol and 30 % distilled water
- · Soft, lint-free cloth

Prerequisites

- The device is switched off and disconnected from the mains/power line.
- The device is defrosted.
- 1. Moisten the lint-free cloth with decontamination agent.
- 2. Clean the surfaces using a lint-free cloth.

The surfaces are coated with a decontamination agent.

- 3. Allow the decontamination agent to take effect.
- 4. Wipe off the decontamination agent with deionized water.
- 5. Allow the surfaces to dry.

7.4 Fuses

Fuses may only be replaced by authorized service technicians. Users must not replace the fuses.

7.5 Checking the alarm

7.5.1 Checking indicator lights and the signal tone

▶ Press the alarm test/mute softkey.

As long as the **alarm test/mute** softkey is pressed, all indicator lights are lit. The signal tone sounds. The display shows 8888.

7.5.2 Checking the alarm in the case of a mains/power outage

Prerequisites

- The back-up circuit is activated.
- ▶ Switch off the device using the mains/power switch.

The **power fail** indicator light lights up.

The display shows the interior temperature and flashes at intervals of 10 s.

The signal tone sounds on the device.

If the device is connected to a building management system via the remote alarm interface, the alarm is forwarded to the building management system.

7.6 Safety checklist

- 1. Fill in the safety checklist prior to repair or service of the device.
- 2. Hand a copy of the safety checklist to the authorized service technician.

				eppendor
1.	Freezer contents	□Yes	□No	
	Risk of infection	□Yes	□No	
	Risk of toxicity	□Yes	□No	
	Risk from radioactive sources	□Yes	□No	
	(List all potentially hazardous materials that have Notes:	ave been s	stored in this unit.)	
2	Contamination of the unit:			
-	Unit interior	□Yes	□No	
	No contamination	□Yes		
	Decontaminated	□Yes		
	Contaminated	□Yes		
	Others			
3.	Instructions for safe repair/maintenance of the a) The unit is safe to work on b) There is some danger (see below) Procedure to be adhered to in order to reduce	□Yes □Yes	□No	w.
Sig Add	te : gnature : dress, Division : lephone :			
Pro	oduct name : odel : rial number :			
Мо				
Mo Sei	te of installation :			
Mo Sei Dai	te of installation : ease decontaminate the unit yourself before cal	lling the se	rvice engineer.	

8 Troubleshooting

8.1 General errors

If you are unable to resolve the error with the suggested measures, please contact your Eppendorf partner. The address can be found on our website: www.eppendorf.com.

8.1.1 Outer door

Problem	Cause	Solution
The outer door cannot be opened.	The door handle is locked.	▶ Unlock the door handle.
	The <i>auto vent</i> valve is blocked. The negative pressure in the interior is preventing the outer door from being opened.	 Wait until pressure compensation has taken place. Balancing the pressure takes 1 to 2 hours. After opening the outer door, remove the ice from the <i>auto vent</i> valve.
The keyboard is not responding.	The keyboard is faulty.	Switch the device off and back on.Contact your Eppendorf partner.

8.1.2 Air filter

Problem	Cause	Solution
The filter clean indicator light lights up.	The air filter is contaminated.	▶ Clean the air filter.

8.2 Software error messages

Problem	Cause	Solution
Code <i>E-01</i> The device triggers an alarm.	The temperature sensor for measuring the interior temperature does not work.	► Contact your Eppendorf partner.
Code <i>E-02</i> The device triggers an alarm.	The temperature sensor at the condenser does not work.	► Contact your Eppendorf partner.
Code <i>E-03</i> The device triggers an alarm.	The temperature sensor at the heat exchanger does not work.	► Contact your Eppendorf partner.
Code <i>E-04</i>	The temperature at the condenser is too high.	 Clean the air filter. Establish an ambient temperature that complies with the technical data. Contact your Eppendorf partner.

Problem	Cause	Solution
Code <i>E-05</i>	The temperature sensor for measuring the ambient temperature does not work.	► Contact your Eppendorf partner.
Code <i>E-06</i> – <i>E12</i> , <i>E14</i> – <i>E20</i>	Internal error messages	► Contact your Eppendorf partner.
Code E-13	The fan does not work.	► Contact your Eppendorf partner.

8.3 Mains/power outage

During a mains/power outage, no mains/power supply is supplied to the device. The device will trigger the "Mains/power outage" alarm . All danger signals go out once the device is supplied with power again.

The interior temperature may rise during a longer mains/power outage.

- If the interior temperature is below the alarm limit for the maximum temperature in the interior after the mains/power outage, the device continues to operate normally.
- If the interior temperature is above the alarm limit for the maximum interior temperature, the "Interior temperature" alarm is triggered after the delay time has elapsed.

8.4 Heating up of the interior

In case of a mechanical or electrical defect, the temperature inside the device may rise after some time. The temperature inside the device rises when the doors of the device are open and warm ambient air enters the device.

When the temperature inside the device exceeds the alarm limit, the "Interior temperature" alarm is triggered.

Open the outer doors and inner doors as briefly as possible to avoid that the temperature inside the device rises. In case of a malfunction, the back-up systems keeps the interior temperature stable over a longer period of time.

9 Transport, storage and disposal

9.1 Decommissioning

Tools and auxiliary equipment

· Adhesive tape

Prerequisites

- Racks and samples have been transferred to another ULT freezer.
- 1. Secure the air intake grille with adhesive tape.
- 2. Disable the back-up circuit (see p. 40).
- 3. Disconnect the device from the voltage supply (see p. 40).
- 4. Remove the safety clamp of the mains/power cord. Remove the mains/power cord from the device.
- 5. Defrost the device (see p. 49).
- 6. Decontaminate the device (see p. 52).

9.2 Transport



DANGER! Risk of severe injury from tipping the device over during transport

If the device tips over and falls on someone, that person sustains fatal injuries.

- ▶ Transport the device with a sufficient number of helpers.
- ▶ Observe the transport instructions in the operating manual.



CAUTION! Risk of foot injury due to little ground clearance

Feet can easily get trapped under the device.

Wear safety shoes with steel toes.



CAUTION! Risk of cut injuries due to sharp-edged components

When working on the air intake grille and on the leveling feet, cut injuries can be caused by sharp-edged components on the bottom of the device.

▶ Wear protective gloves to avoid cut injuries.



NOTICE! Risk of device damage due to lifting the device without the original pallet

Lifting the device without the original pallet will damage the base of the device.

- 1. Place the device on the original pallet.
- 2. Secure the device.
- 3. Use a transport aid to lift the device.



NOTICE! Damage to the compressor and refrigeration cycle during transport

Tilting the device or transporting it in a horizontal position will damage the compressors and the refrigeration cycle. Refrigerant and oil may leak out.

Shocks may dislodge the compressors from the brackets.

▶ Transport the device in upright position.

- ▶ Move the device with due caution and care. Do not knock the device into anything.
- ▶ Protect the device from impacts.
- ▶ After setting up the device, wait for 6 h before switching it on.



NOTICE! Damage to the door handle due to too high loads

Pulling or pushing the device on the door handle during transport may damage the door handle.

• Grip the device at the housing to pull or push it.



NOTICE! Transport damage to the control panel

The control panel protrudes from the door. The control panel may become damaged during transport.

▶ Only transport the device with the transport protection attached to the control panel.

9.2.1 Preparing the device for transport

Personal protective equipment

· Protective gloves

Tools and auxiliary equipment

- · Open-end wrench
- · Adhesive tape

Prerequisites

- Racks and samples have been transferred to another ULT freezer.
- 1. Secure the air intake grille with adhesive tape.
- 2. Attach the transport protection to the control panel.
- 3. Disable the back-up circuit (see p. 40).
- 4. Disconnect the device from the voltage supply (see p. 40).
- 5. Remove the safety clamp of the mains/power cord. Remove the mains/power cord from the device.
- 6. Use the open-end wrench to turn the leveling feet upward (Fig. 4-1 on p. 31).

9.2.2 Transporting the device

Personal protective equipment

• Protective clothing, safety shoes

Tools and auxiliary equipment

- Transport aid
- Original pallet

General transport

- 1. Transport the device in an upright position.
- 2. Grasp the device by the housing and wheel it to its new location. Do not grasp the device by the door handle.

Lifting the device

- 3. Place the device on the original pallet and secure it.
- 4. Lift the device with a transport aid.

Sloping surfaces

- 5. Transport the device sideways over ramps.
- 6. Do not transport the device over ramps at an angle $> 17 \% (10^{\circ})$.

Narrow passageways

- 7. Open the door of the device 180°.
- 8. Push the device, with one side panel to the front, through the passageway.

 If the passageway is too narrow for the device, housing parts may have to be disassembled. Contact your local Eppendorf partner for more information.

Outside of buildings

9. Transport the device with a transport aid.

The heavy-duty castors are not suitable for transport outside of buildings.



If you need help to transport the device, contact the authorized service.

9.3 Shipment

9.3.1 Shipping regulations

ULT freezers that contain more than 100 g of flammable refrigerant are classified as refrigerating machines containing flammable, non-toxic, liquefied gas (UN no. 3358).

The ULT freezer contains more than 100 g of flammable refrigerant and must not be transported via air freight.

9.3.2 Shipping the device



WARNING! Risk of personal injury due to contamination.

People may get contaminated if you store or ship a contaminated device.

▶ Clean and decontaminate the device before shipping or storage.



NOTICE! Risk of damage due to incorrect packing.

Eppendorf AG is not liable for any damage caused by improper packing.

- ▶ Only store and transport the device in its original packing.
- If you do not have the original packing, request the original packing from Eppendorf AG.

Prerequisites

- The device has been taken out of operation.
- The device has been cleaned and decontaminated.
- The original packing is available.
- 1. Download the "Decontamination declaration for product returns" from the www.eppendorf.com website.
- 2. Complete the decontamination certificate.
- 3. Pack the device.
- 4. Put the decontamination certificate into the packing.
- 5. Ship the device according to the shipment regulations.



For service and repairs, send the device to Eppendorf AG or to an authorized service partner.

9.4 Disposal

The product must be disposed of in accordance with the relevant legal provisions.

Information on the disposal of electrical and electronic devices in the European Community:

Within the European Community, the disposal of electrical devices is regulated by national regulations based on EU Directive 2012/19/EU pertaining to waste electrical and electronic equipment (WEEE).

According to these regulations, any devices supplied after August 13, 2005, in the business-to-business sphere, to which this product is assigned, may no longer be disposed of in municipal or domestic waste. They are marked with the following symbol to indicate this:



Do not dispose of batteries as household waste. Dispose of batteries in accordance with the local regulations.

As the disposal regulations may differ from one country to another within the EU, please contact your supplier for more information.

Transport, storage and disposal CryoCube® F740h English (EN)

62

10 Technical data

10.1 Power supply

Mains/power supply voltage	100 V - 230 V ±10 %
Mains/power frequency	50 Hz - 60 Hz
Current consumption	100 V – 230 V (50 Hz); 15 A – 6 A 100 V – 220 V (60 Hz); 12 A – 6 A
Electromagnetic compatibility (EMC)	The device meets the following requirements: • IEC/EN 61326-1 • EN 55011 (CISPR 11) • FCC Part 15 – Class A
Overvoltage category	II
Degree of pollution	2


10.2 Ambient conditions

10.2.1 Operation

Ambience	For indoor use only
Ambient temperature	15 °C – 32 °C
Relative humidity	Maximum 80 %, non-condensing
Atmospheric pressure	80 kPa – 106 kPa

10.3 Dimensions

10.3.1 External dimensions

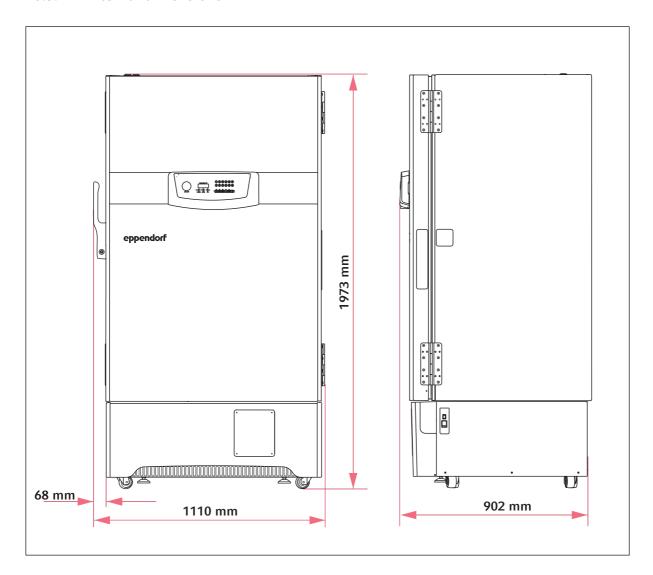

10.3.2 Internal dimensions

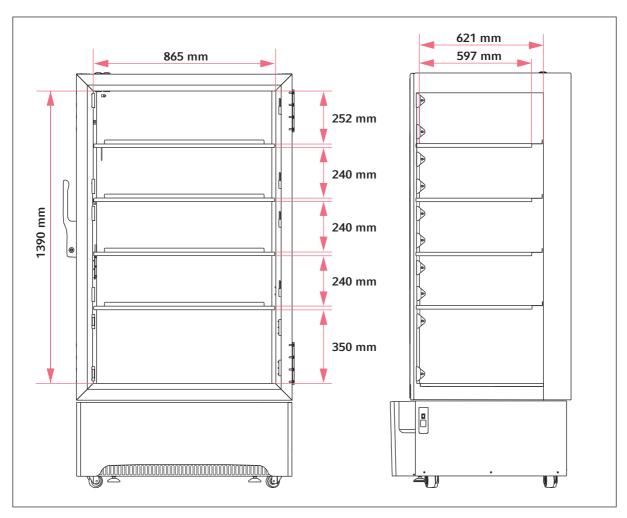


Fig. 10-1: Internal dimensions for devices with 5 inner compartments

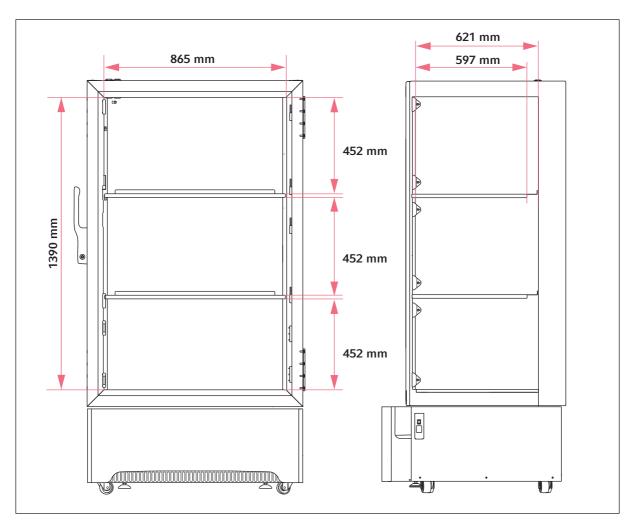


Fig. 10-2: Internal dimensions for devices with 3 inner compartments

10.3.3 Packing dimensions

Width	1200 mm
Depth	1045 mm
Height	2225 mm

10.4 Weight

Device	Model with 3 compartments	Model with 5 compartments
CryoCube F740h	326 kg	342 kg
Packaged device	Model with 3 compartments	Model with 5 compartments
CryoCube F740h	376 kg	392 kg

10.5 Noise level

CryoCube F740h	47.8 dB (A)
----------------	-------------

10.6 Interfaces

BMS (remote alarm)	24 V, 1 A
Serial interface	RS-485 (for internal use)

Only connect devices that comply with the IEC 60950 (UL 60590) standards to the interfaces.

10.7 Temperature control

10.7.1 Temperature range

Setting range	-50 °C to -86 °C
---------------	------------------

10.7.2 Times for cooling and heating the interior

Cooling from 20 °C to -80 °C	CryoCube F740h (230 V)	3 h 45 min
Heating from -80 °C to 0 °C The device is 2/3 full.	CryoCube F740h (230 V)	41 h

10.7.3 Cooling of the refrigeration cycle

CryoCube F740h	Air cooling

10.7.4 Refrigerant

Device	Refrigeration cycle 1	Refrigeration cycle 2
CryoCube F740h	R-290 (96 g)	R-170 (106 g)

Refrigerant	Components
R-170	Ethane
R-290	Propane

10.8 Additional specifications

10.8.1 Capacity and carrying capacity

Capacity	740 L
Carrying capacity per inner shelf	150 kg (with an evenly distributed load)
Carrying capacity per device	420 kg

10.8.2 Materials

Outer door insulation	Vacuum insulation panels Polyurethane foam
Device insulation	Vacuum insulation panels Polyurethane foam
Interior	Stainless steel (304 2B)

11 Ordering information

11.1 Accessories

11.1.1 Back-up systems

Order no.	Description
(International)	
	CO2 back-up system
F652999005	100 V/50 Hz - 60 Hz
U9043-0002	120 V - 220 V/60 Hz
U9043-0004	230 V/50 Hz
	LN2 back-up system
F652999006	100 V/50 Hz - 60 Hz
U9044-0002	120 V - 220 V/60 Hz
U9044-0004	230 V/ 50 Hz

11.1.2 Chart recorder

Order no.	Description
(International)	
	Chart recorder type 2
	Connection to mains/power supply in the ULT freezer
F652999001	100 V/120 V, 50 Hz – 60 Hz
F652999002	208 V – 230 V, 50 Hz – 60 Hz
	Discs for chart recorder type 2
	-100 °C - 0 °C
F652999003	60 pieces
	Pens for chart recorder type 2
F652999004	3 pieces

11.1.3 Racks for devices with 3 compartments

Order no.	Description
(International)	
	Rack with drawers
	material stainless steel
6001 072.210	for 32 boxes, box height 53 mm
6001 072.910	for 24 boxes, box height 63.5 mm
6001 072.310	for 20 boxes, box height 78 mm
6001 072.410	for 16 boxes, box height 103 mm
6001 072.510	for 12 boxes, box height 133 mm
	Rack with side access
	material stainless steel
6001 071.210	for 32 boxes, box height 53 mm
6001 071.910	for 24 boxes, box height 63.5 mm
6001 071.310	for 20 boxes, box height 78 mm
6001 071.410	for 16 boxes, box height 103 mm
6001 071.510	for 12 boxes, box height 133 mm
6001 071.110	with compartments for 48 deepwell plates up to 53 mm

11.1.4 Racks for the compartments 1 – 4 for devices with 5 compartments

Order no.	Description			
(International)				
	Rack with drawers			
	material stainless steel			
6001 022.210	for 16 boxes, box height 53 mm			
6001 022.910	for 12 boxes, box height 63 mm			
6001 022.310	for 8 boxes, box height 78 mm			
6001 022.410	for 8 boxes, box height 103 mm			
	Rack with side access			
	material stainless steel			
6001 021.210	for 16 boxes, box height 53 mm			
6001 021.910	for 12 boxes, box height 63 mm			
6001 021.310	for 8 boxes, box height 78 mm			
6001 021.410	for 8 boxes, box height 103 mm			
6001 021.110	with compartments for 24 deepwell plates up to 53 mm			

11.1.5 Racks for compartment 5 for devices with 5 compartments

Order no.	Description			
(International)				
	Rack with drawers			
	material stainless steel			
6001 082.210	for 24 boxes, box height 53 mm			
6001 082.910	for 20 boxes, box height 63.5 mm			
6001 082.310	for 16 boxes, box height 78 mm			
6001 082.410	for 12 boxes, box height 103 mm			
6001 082.510	for 8 boxes, box height 133 mm			
	Rack with side access			
	material stainless steel			
6001 081.210	for 24 boxes, box height 53 mm			
6001 081.910	for 20 boxes, box height 63.5 mm			
6001 081.310	for 16 boxes, box height 78 mm			
6001 081.410	for 12 boxes, box height 103 mm			
6001 081.510	for 8 boxes, box height 133 mm			
6001 081.110	with compartments for 36 deepwell plates up to 53 mm			

11.1.6 Cardboard boxes and box dividers

Order no.	Description			
(International)				
	Cardboard box			
B50-SQ	width 133 mm, depth 133 mm, height 50 mm			
B75-SQ	width 133 mm, depth 133 mm, height 75 mm			
B95-SQ	width 133 mm, depth 133 mm, height 100 mm			
	Box divider			
D49	for 7 × 7 vessels, maximum vessel diameter 17.4 mm			
D64	for 8 × 8 vessels, maximum vessel diameter 15 mm			
D81	for 9 × 9 vessels, maximum vessel diameter 13 mm			
D100	for 10 × 10 vessels, maximum vessel diameter 11.8 mm			

11.1.7 Eppendorf Storage Boxes

Order no.	Description
(International)	
	Eppendorf Storage Box 10 × 10, 2 inch
	height 52.8 mm, for 100 cryogenic tubes with internal thread
0030 140.508	3 pieces
	Eppendorf Storage Box 9 x 9, 2 inch
	height 52.8 mm, for 81 screw cap (cryogenic) tubes 1 mL - 2 mL
0030 140.516	3 pieces
	Eppendorf Storage Box 8 × 8, 2 inch
	height 52.8 mm, for 64 micro test tubes 1 mL - 2 mL
0030 140.524	3 pieces
	Eppendorf Storage Box 5 × 5, 2.5 inch
	height 63.5 mm, for 25 micro test tubes 5 mL
0030 140.532	4 pieces
	Eppendorf Storage Box 9 x 9, 3 inch
	height 76.2 mm, for 81 screw cap (cryogenic) tubes 3 mL
0030 140.540	2 pieces
	Eppendorf Storage Box 9 x 9, 4 inch
	height 101.6 mm, for 81 screw cap (cryogenic) tubes 4 mL - 5 mL
0030 140.567	2 pieces
	Eppendorf Storage Box 5 x 5, 5 inch
	height 127 mm, for 25 conical tubes 15 mL
0030 140.583	2 pieces
	Eppendorf Storage Box 3 x 3, 5 inch
	height 127 mm, for 9 conical tubes 50 mL and 4 conical tubes 15 mL
0030 140.591	2 pieces
	Eppendorf Storage Box 5 x 5, 3 inch
	height 76.2 mm, for 25 screw cap tubes 5 mL
0030 140.613	2 pieces

Index	Operation63
	Technical data
Α	Mains/power supply63
Alarm24	W
В	Weight66
Battery switch Switching off	
D	
Degree of pollution63	
Device Switching off	
Disposal61	
E	
Electromagnetic compatibility (EMC)63	
Electionagnetic compatibility (EMC)	
М	
Mains/power supply63	
N	
Noise level67	
Noise level	
s	
Switching off Back-up circuit	
Switching on Back-up circuit	
Т	
Technical data	
Ambient conditions63	
Noise level 67	

Index CryoCube® F740h English (EN)

74

eppendorf

Declaration of Conformity

The product named below fulfills the requirements of directives and standards listed. In the case of unauthorized modifications to the product or an unintended use this declaration becomes invalid. This declaration of conformity is issued under the sole responsibility of the manufacturer.

Product name:

CryoCube® F740h, CryoCube® F740hi, CryoCube® F740hiw

including accessories

F740320011 F740320021 F740320031 F740320041 F740320111 F740320131 F740340011 F740340021 F740340041

Product type:

Ultra-low temperature freezer

"i" designates model with touch user interface, in general
"h" designates model with refrigerant hydrocarbon, air-cooled
"hi" designates model with refrigerant hydrocarbon, air-cooled
"hiw" designates model with refrigerant hydrocarbon, water-cooled

Relevant directives / standards:

2014/35/EU: EN 61010-1, IEC 61010-1

UL 61010A-1, CAN/CSA C22.2 No. 61010-1

2014/30/EU: EN 61326-1

47 CFR FCC part 15

2011/65/EU: EN 50581

2006/42/EC: EN 378-2 (partial)

Further applied

standards: UL 60730-1, UL 471, CAN/CSA-E60730-1, CSA C22.2 No. 120,

IEC 60335-2-89

Hamburg, October 15, 2018

Dr. Wilhelm Plüster Management Board

Dr. Sven Bülow Head of Business Unit Sample Management

Your local distributor: www.eppendorf.com/contact Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany eppendorf@eppendorf.com ISO 9001 Certified ISO 13485 Certified ISO 14001 Certified

CERTIFICATE OF COMPLIANCE

Certificate Number 2018-03-29; 2018-07-09 (A1)-E215059

Report Reference E215059-D1001-1/A1/C0-UL-UL

Issue Date 2018-03-29; 2018-07-09 (A1)

Issued to: Eppendorf A G

Applicant Company: Barkhausenweg 1

22339 Hamburg Germany

Listed Company: Same as Applicant

This is to certify that Laboratory Freezer

representative samples of CryoCube F740, CryoCube F740i, CryoCube F740iw, CryoCube

F740h, CryoCube F740hi, CryoCube F740hiw

Have been investigated by UL in accordance with the

Standard(s) indicated on this Certificate.

Standard(s) for Safety: UL 61010-1, 3rd Edition, May 11, 2012, Revised July 15 2015,

CAN/CSA-C22.2 No. 61010-1-12, 3rd Edition, Revision dated

July 2015

Additional Standards: UL 471 (edition Tenth 06/17/2016), Sections 42, 44, 49 and 68

and CSA C22.2 No. 120-13 (March 2013) Sections 6.11, 6.3, 6.4

and 6.26

Models F740hi, F740hiw and F740h were investigated to UL 471 Tenth Edition, with revisions through December 8, 2016, and the Canadian Standard for Refrigeration Equipment, CAN/CSA C22.2 No. 120-13 dated March 2013 under report SA45018

Vol1, Sec. 1.

Fluid-containing parts of refrigeration systems meet the relevant

pressure-related requirements of IEC 60335-2-89.

Additional Information: See the UL Online Certifications Directory at

www.ul.com/database for additional information.

Helena y. wol

Only those products bearing the UL Certification Mark should be considered as being covered by UL's Certification and Follow-Up Service.

Look for the UL Certification Mark on the product.

This is to certify that representative samples of the product as specified on this certificate were tested according to the current UL requirements.

ruce Mahrenholz, Assistant Chief Engineer, Global Inspection and Field Services, UL LLC
Helena Y. Wolf, Director, Global Market Access Operations, UL LLC
pseph Hosey, General Manager, Director of Sales – Canada, UNDERWRITERS LABORATORIES OF CANADA INC.

Any information and documentation involving UL Mark services are provided on behalf of UL LLC (UL) or any authorized licensee of UL. For questions, please contact a local UC Customer Service Representative www.ul.com/contactus



Evaluate Your Manual

Give us your feedback. www.eppendorf.com/manualfeedback